

Moisture Determination of Cement Raw Mix

Precisa moisture analysis XM60/EM120 Halogen



Moisture Content



160°C



Sample Weighing



Auto-Stop



This information is intended as a guideline for development of a moisture analysis method and parameter setting.

Materials vary throughout the world and we will be more than happy to test customers' individual samples.

Description:	Cement raw mix, fine powder
Utensils:	Aluminium pan
Preparation:	Spread thinly on the pan
Sample weight (+/- 10%):	2g
Temperature program:	Standard
Temperature:	105 °C
Standby Temperature	100 °C
Stop-Criteria:	Autostop 2d / 60 s
Moisture content: (average of 4 measurements)	0.68%
Standard Deviation (+/-):	0.04%
Average analysis time:	1.8 min